


Aspergus medium for Cryptomyces

(1) Broth :-

- A { Peptone (Bacto-peptone) 10.
Glucose 20.
Distilled water about 300. cc

Divide in water bath in suitable large flask

- B { Description of aspergus -
Take a tin of aspergus heads
Boil in a pan with about 500 cc distilled water till
nicely extracted
Filter through cheesecloth
The fairly clear solution is then added to A

- C - Adjust pH to 7.0 (or 7.2) (old method of titration)
Add 50 cc glycerin
& make up to one liter
Autoclave & keep in 100 cc flasks: 
(or less)

Use one for the ^{stock} subculture kept covered growth

(2) Agar :-

Add to a given volume of (1) to give 2-3 % agar
For slope tubes & for stab tubes (used for plate pouring).